
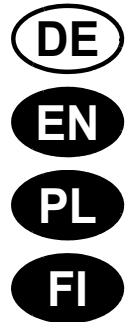
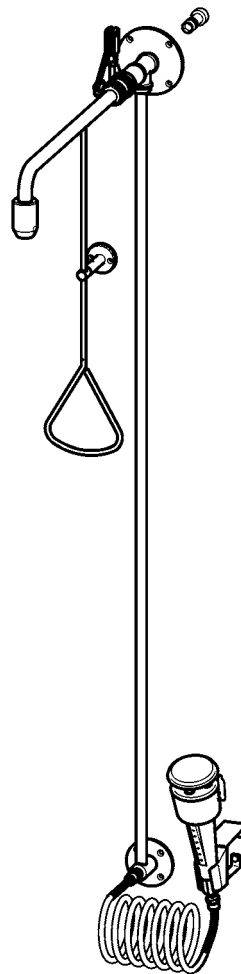
Best.-Nr.: 2030019212  
FAID0010

Best.-Nr.: 2030019221  
FAID0011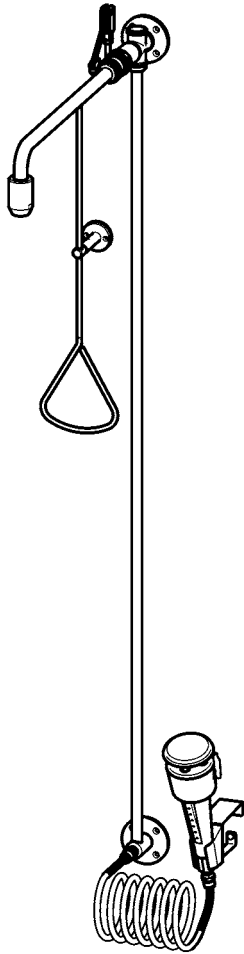

<b>EN</b>	.....	<b>3</b>
-----------	-------	----------

Please refer to the graphics in the German Installation and Operating Instructions.

<b>PL</b>	.....	<b>12</b>
-----------	-------	-----------

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Emergency shower combination

Please refer to the graphics in the German installation and operating instructions.

## Table of contents

1. Abbreviations and units . . . . .	3
2. Key . . . . .	4
3. Warranty . . . . .	4
4. Important notes . . . . .	4
<b>Description of product</b>	
5. Application . . . . .	4
6. Technical specifications . . . . .	5
7. Scope of delivery . . . . .	6
8. Dimensions . . . . .	6
9. Installation example . . . . .	6
<b>Assembly, function and commissioning</b>	
10. Installation of body shower water connection from above . . . . .	6
11. Installation of body shower water connection from behind . . . . .	7
12. Installation of eye shower . . . . .	9
13. Function . . . . .	9
<b>Maintenance</b>	
14. Maintenance and care . . . . .	10
15. Replacing a filter . . . . .	10
16. Replacing the flow rate regulator . . . . .	10
17. Replacing the non-return valve . . . . .	10
18. Troubleshooting . . . . .	11
19. Spare parts . . . . .	11

## 1. Abbreviations and units

Best.-Nr.	Aquarotter order number
Conversion	1 mm = 0.03937 inches 1 inch = 25.4 mm

All length specifications in the graphics are in mm.

## 2. Key

---

### **Warning!**

Failure to observe can result in bodily injury or even death.

### **Caution!**

Failure to observe can result in material damage.

### **Important!**

Failure to observe can cause the product to malfunction.

Useful information for optimum handling of the product.

## 3. Warranty

---

Liability is accepted in accordance with the General Terms and Conditions of Business and Supply.

Use only genuine spare parts!

## 4. Important notes

---

- Installation, commissioning and maintenance must be performed only by a qualified expert in accordance with the instructions provided, legal requirements and recognised engineering standards.
- All technical connection regulations specified by the local water and electricity supply companies must be observed.
- We reserve the right to make changes.

## 5. Application

---

Emergency showers are prescribed first aid installations for work stations where hazardous materials are being handled.

## 6. Technical specifications

---

### Body shower

Minimum flow pressure: 0.5 bar

Maximum stagnation pressure: 10 bar

Volumetric flow:

0.5 bar flow pressure: 45 l/m

1.0 bar flow pressure: 65 l/m

1.5 bar flow pressure: 80 l/m

2.0 bar flow pressure: 90 l/m

2.5 bar flow pressure: 100 l/m

3.0 bar flow pressure: 110 l/m

Assembly height: Lower edge of shower head ( $220 \pm 10$ ) cm above floor as per EN 15154 Part 1

### Eye shower

Version: Hand shower plastic green/white  
Wall mounting, tilted 20° forwards, spiral hose and protective cap with user information

Minimum flow pressure: 1 bar

Flow pressure range: 1-4 bar  
This range meets the requirements of EN 15154.

Volumetric flow: Constant 6-9 l/s at a flow pressure of 1 bar

## 7. Scope of delivery

Pos.	Name	FAID0010	FAID0011
1	Shower head	X	X
2	Wall bracket	X	X
3	Ball valve	X	X
4	Hand actuator and guide post support with plugging option	X	X
5	Connecting pipe	X	X
6	Wall plate with T-piece, reducer	X	X
7	Wall plate with cross piece	X	
8	Wall connection and plug socket		X
9	Accessories kit: First aid sign, Mounting set	X	X
10	Eye shower with safety cap	X	X
11	Spiral supply hose with filter	X	X
12	Wall mounting with attachment set	X	X
13	First aid sign	X	X
	Installation and operating instructions	X	X

## 8. Dimensions

For special designs, see the customer drawing for dimensions

## 9. Installation example

## 10. Installation of body shower water connection from above

### Important!

- Prior to the assembly process, flush the pipework in accordance with DIN 1988.
- Mount the emergency shower in accordance with EN 15154 Part 1.

 The manual actuator of the emergency shower can be plugged.

 To protect the tiles, use masking tape when marking and drilling. Drill at low speed.

 Depending on the composition of the wall, special wall plugs may have to be used (to be provided by the customer).

**10.1** Screw the connecting pipe (b) into the wall plates so that a seal is formed.

a: Wall plate with cross piece

c: wall plate with T-piece

### **Important!**

Pay attention to precisely vertical alignment during installation.

- 10.2 Align the wall panel (a)  $2310 \pm 100$  mm above the finished floor in accordance with the water intake, and mark the drill holes.
- 10.3 Drill in accordance with these markings.
- 10.4 Insert the wall plugs.
- 10.5 Screw on the wall plates.  
a: Wall plate with cross piece  
c: wall plate with T-piece
- 10.6 Screw in the closed ball valve as far as possible, creating a seal and bringing it into position as follows:
  - The valve axis must point to the left.
  - The pivot motion of the valve lever must be aligned so that it is precisely vertical.
- 10.7 Secure the ball valve (d) with the two grub screws (c).
- 10.8 Place the wall support (e) onto the guide (g) on the manual actuator and secure with the grub screw (f).
- 10.9 Place the valve lever into the "close" position.  
A: Closed  
B: Open
- 10.10 Screw the manual actuator (j) into the fork head (h) of the deflection lever and lock with a nut.
- 10.11 Using the spirit level, align the manual actuator (j) so that it is vertical.
- 10.12 Mark drill holes for the wall support (e) 40 mm above the triangular handle.
- 10.13 Drill in accordance with these markings.
- 10.14 Insert the wall plugs.
- 10.15 Screw on the wall support (e).
- 10.16 Mount the wall arm in a sealing manner and bring it into position.
- 10.17 Screw the shower head (l) onto the wall arm (k) so that it seals.

## **11. Installation of body shower water connection from behind**

---

### **Important!**

- Prior to the assembly process, flush the pipework in accordance with DIN 1988.
- Mount the emergency shower in accordance with EN 15154 Part 1.

-  The manual actuator of the emergency shower can be plugged.
-  To protect the tiles, use masking tape when marking and drilling. Drill at low speed.
-  Depending on the composition of the wall, special wall plugs may have to be used (to be provided by the customer).

- 11.1 Screw the connecting pipe (b) into the wall plates so that a seal is formed.  
a: Wall plate with cross piece  
c: wall plate with T-piece
-  **Important!**  
The internal thread G  $\frac{3}{4}$  for the locally provided water connection may not protrude from the wall by more than 1 mm, and may not extend into the wall by more than 15 mm.
- 11.2 Place the water connection centrally at a height of  $2310 \pm 100$  mm above the finished floor.
- 11.3 Screw in the plug-in socket ( ) with an Allen key to form a seal.
-  **Important!**  
Pay attention to precisely vertical alignment during installation.
- 11.4 Slide the wall plate (a) onto the plug-in socket (c) and align it.
- 11.5 Mark the drill holes.
- 11.6 Drill in accordance with these markings.
- 11.7 Insert the wall plugs.
- 11.8 Slide the wall plate (a) onto the plug-in socket (c) as far as the wall and fix in place with the grub screw (d).
- 11.9 Screw on the wall plates.  
a: Wall plate with cross piece  
c: wall plate with T-piece
- 11.10 Screw in the closed ball valve as far as possible, creating a seal and bringing it into position as follows:
- The valve axis must point to the left.
  - The pivot motion of the valve lever must be aligned so that it is precisely vertical.
- 11.11 Secure the ball valve (f) with the two grub screws (e).
- 11.12 Place the wall support (g) onto the guide (j) on the manual actuator and secure with the grub screw (h).
- 11.13 Place the valve lever into the "close" position.  
A: Closed  
B: Open
- 11.14 Screw the manual actuator (l) into the fork head (h) of the deflection lever and lock with a nut.
- 11.15 Using the spirit level, align the manual actuator (l) so that it is vertical.
- 11.16 Mark drill holes for the wall support (g) 40 mm above the triangular handle.
- 11.17 Drill in accordance with these markings.
- 11.18 Insert the wall plugs.
- 11.19 Screw on the wall support (g).
- 11.20 Mount the wall arm in a sealing manner and bring it into position.
- 11.21 Screw the shower head (n) onto the wall arm (m) so that it seals.

## 12. Installation of eye shower

---

### **Important!**

- Prior to the assembly process, flush the pipework in accordance with DIN 1988.
- Mount the eye shower in accordance with EN 15154 Part 2.

12.1 Mount the wall bracket (b) to the extension (a).

12.2 Mark the drill holes for the wall bracket at a height (c) of 590 mm to 990 mm from the upper edge of the finished floor.

12.3 Drill holes in accordance with these markings.

12.4 Insert the wall plugs.

12.5 Mount the extension (a).

### **Caution!**

Note the installation position of the filter.

12.6 Insert the filter (b) into the hose (c).

12.7 Connect the eye shower (a) to the hose (c).

12.8 Insert the hose (c) into the wall bracket (e).

12.9 The handle bar (d) must face forwards.

12.10 Screw the reducer (g) into the wall plate (f) to seal it.

12.11 Screw the hose (c) with the seal onto the reducer (g).

12.12 Establish the water connection.

12.13 Check the water connection for leaks.

12.14 Plug the manual actuator if need be.

## 13. Function

---

13.1 Pull the manual actuator.

- The water will flow.

 The valve is not self-closing!

13.2 Place the manual actuator back in its start position.

- The water stops flowing.

### **Stationary use**

13.3 Press the trigger button.

- The water will flow.
- The safety cap will fall off.

13.4 Press the red stop button.

- The water flow will stop.

 The valve is not self-closing! The user can use both hands to keep their eyes open!

### **Mobile use**

 Depending on the length of the hose, the eye shower can also be pulled out of its bracket.

## 14. Maintenance and care

---

Check the emergency shower combination once a month to make sure that it is in working order. (Guidelines for Laboratories BGI/GUV-I 850-0)

As a measure for the prevention of microbiological contamination, it is recommended that you replace the fill water in the fitting at shorter intervals by opening the valve.

Use suitable cleaning agents which do not attack the tap in a proper manner and rinse with water after use. High-pressure cleaners must not be used for cleaning.

## 15. Replacing a filter

---

15.1 Close the water supply.

15.2 Take the eye shower (a) out of the wall bracket.

15.3 Loosen the eye shower (a) from the hose (d).

15.4 Replace the (c) filter.

### **Important!**

Note the installation position of the filter.

The handle bar (b) must face forwards.

15.5 Reassemble in reverse order.

## 16. Replacing the flow rate regulator

---

16.1 Interrupt the water supply.

16.2 Take the eye shower (a) out of the wall bracket.

16.3 Loosen the hose (b).

16.4 Use a bent wire to lift out the flow volume regulator and replace it.

### **Caution!**

Note the installation position of the non-return valves.

Do not forget the seal.

16.5 Reassemble in reverse order.

## 17. Replacing the non-return valve

---

### **Caution!**

The non-return valve must be replaced every 2 years.

17.1 Loosen the screw (a).

17.2 Remove the shower base (b).

17.3 Remove the spray heads (c) and the seals (d).

17.4 Unscrew the attachment part (e).

17.5 Remove the disc (f).

17.6 Push out the non-return valve (g).

17.7 Reassemble in reverse order.

## 18. Troubleshooting

Fault	Cause	Remedy
Water does not flow	– Water supply interrupted	⇒ Restore it
Manual actuator is catching	– Ball valve not correctly aligned	⇒ Align it
	– Pull rod or manual actuator bent	⇒ Straighten it
	– Ball valve stiff	⇒ Replace it
Body shower is dripping	– Ball valve defective	⇒ Replace it
Water volume too low	– Flow volume regulator faulty	⇒ Replace it
	– Dirty filter	⇒ Clean it

If you cannot rectify a fault or if it is not in the list, please contact our customer service department!

## 19. Spare parts

Description	Order No.	
1 Hand shower complete . . . . .	2000105863	7 Filter (packaging unit 10x) . . . . . 2000104732
2 Spray head (packaging unit 4x) . . . . .	2000104725	Service kit consisting of:
3 Safety cap with ball chain . . . . .	2000105849	O-rings . . . . . 2000105864
4 Non-return valve (packaging unit 10x) . . . . .	2000104744	Ball valve . . . . . 2030025022
5 Spiral connection hose . . . . .	2000105844	
6 Flow volume regulator . . . . .	2000105842	

 Other spare parts on request; see the wall bracket for material numbers.

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KWC ME LLC Ras Al Khaimah,
KWC DVS Ltd - Northern Office
KWC DVS Ltd - Southern Office
KWC Aquarotter GmbH
KWC Nordics Oy
**KWC Austria GmbH**